

Human Proinsulin ELISA Kit

(Catalog Number: 31420)

For the quantitative determination of human proinsulin
concentrations in serum or plasma

IMD (Hong Kong)

Address: Unit 513, 5/F, Biotech Centre 2, No. 11 Science Park West Avenue,
Hong Kong Science Park, Sha Tin, Hong Kong

Website: www.immunodiagnostics.com.hk

Email: info@immunodiagnostics.com.hk

Tel: (+852) 3502 2780

IMD (Canada)

Address: 3330 Bayview Avenue, Block #6, Toronto, M2M 3R8, Ontario, Canada

Email: info@immunodiagnostics.ca

Tel: +1-437-886-5136

Website: www.immunodiagnostics.com.hk

E-mail: info@immunodiagnostics.com.hk(HK) / info@immunodiagnostics.ca(Canada)

Tel: +852 3502 2780 (HK) / +1-437-886-5136 (Canada)

Version 09/25

TABLE OF CONTENTS

Contents	Page
INTRODUCTION	1
PRINCIPLE OF THE ASSAY	1
INTENDED USE	1
REAGENTS SUPPLIED	1
OTHER MATERIALS REQUIRED, BUT NOT PROVIDED	1
STORAGE	2
PREPARATION OF REAGENTS	2
PREPARATION OF STANDARDS	2
ASSAY PROCEDURE	3
CALCULATION	3
TYPICAL STANDARD CURVE	4
ASSAY CHARACTERISTICS	4
SUMMARY OF ASSAY PROCEDURE	5

Website: www.immunodiagnostics.com.hk

E-mail: info@immunodiagnostics.com.hk(HK) /info@immunodiagnostics.ca(Canada)

Tel: +852 3502 2780 (HK) / +1-437-886-5136 (Canada)

Version 09/25



INTRODUCTION

Proinsulin is a polypeptide of 86 amino acids made in the beta cells of the pancreas and is the precursor molecule for insulin, most proinsulin is converted to insulin and C-peptide secreted into the blood. Increased circulating levels of proinsulin are found in patients with insulinomas, hypoglycemia and hyperinsulinemia.

PRINCIPLE OF THE ASSAY

This assay is a quantitative sandwich enzyme-linked immunosorbent assay (ELISA). The microtiter plate is pre-coated with a monoclonal antibody specific for human proinsulin. Standards, samples and HRP labelled monoclonal antibody specific for human proinsulin are pipetted into the wells and any human proinsulin present is bound by the immobilized antibody. After wash step to remove any unbound reagents, an HRP substrate solution is added and color develops in proportion to the amount of human proinsulin bound initially. The assay is stopped, and the optical density of the wells is determined using a microplate reader. Since the increases in absorbance are directly proportional to the amount of captured human proinsulin, the unknown sample concentration can be interpolated from a reference curve included in each assay.

INTENDED USE

This Human Proinsulin ELISA kit is designed for quantification of human proinsulin in serum and plasma samples.

REAGENTS SUPPLIED

Each kit is sufficient for one 96-well plate and contains the following components:

1. Microtiter Strips (96 wells), coated with a monoclonal antibody against human proinsulin, sealed
2. 10×Wash buffer, 30 mL
3. 1×Assay buffer, 25 mL
4. 100×Detection antibody solution, a HRP labelled monoclonal antibody against human proinsulin, 0.06 mL
5. Human proinsulin standard, 2000 pg of recombinant human proinsulin, lyophilized
6. Substrate solution, 12 mL, ready for use
7. Stop solution, 12 mL, ready for use
8. Plate cover

OTHER MATERIALS REQUIRED, BUT NOT PROVIDED

1. Pipettes and pipette tips
2. 96-well plate or manual strip washer
3. Buffer and reagent reservoirs
4. Paper towels or absorbent paper

Website: www.immunodiagnostics.com.hk

E-mail: info@immunodiagnostics.com.hk(HK) /info@immunodiagnostics.ca(Canada)

Tel: +852 3502 2780 (HK) / +1-437-886-5136 (Canada)



5. Plate reader capable of reading absorbency at 450 nm
6. Distilled water or deionized water

STORAGE

The kit should be stored at 2-8°C upon receipt, and all reagents should be equilibrated to room temperature before use. Remove any unused antibody-coated strips from the human proinsulin microtiter plate, return them to the foil pouch and re-seal. Once opened, the strips may be stored at 2-8°C for up to one month.

PREPARATION OF REAGENTS

Bring all reagents and materials to room temperature before assay.

A. 1×Wash buffer


Prepare 1×Wash buffer by mixing the 10×Wash buffer (30 mL) with 270 mL of distilled water or deionized water. If precipitates are observed in the 10×Wash buffer bottle, warm the bottle in a 37°C water bath until the precipitates disappear. The 1×Wash buffer may be stored at 2-8°C for up to one month.

B. 1×Detection antibody solution

Spin down the 100×Detection antibody solution briefly and dilute the desired amount of the antibody 1:100 with 1×Assay buffer, 50 µL of the 1×Detection antibody solution is required per well. Prepare only as much 1×Detection antibody solution as needed. Return the 100×Detection antibody solution to 2-8°C immediately after the necessary volume is removed.

PREPARATION OF STANDARDS

Human Proinsulin Standards: Reconstitute the lyophilized standard with 1 mL of 1×Assay buffer to generate a standard stock solution of 2000 pg/mL. Allow the standard to sit for 10 minutes with gentle agitation prior to making dilutions. Pipette 500 µL of 1×Assay buffer to 1000, 500, 250, 125, 62.5, 31.2 pg/mL tubes. Use the standard stock solution to produce a serial dilution as shown below.

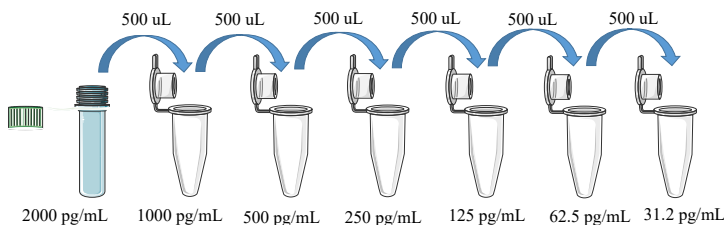

Website: www.immunodiagnostics.com.hk

E-mail: info@immunodiagnostics.com.hk(HK) /info@immunodiagnostics.ca(Canada)

Tel: +852 3502 2780 (HK) / +1-437-886-5136 (Canada)

Version 09/25



1×Assay buffer serves as the zero standard (0 pg/mL). The reconstituted standard stock should be aliquoted and stored at -80°C for one month. Avoid repeating freezing/thawing cycles. Please do not store the diluted standard solutions.

SAMPLE HANDLING

Serum or plasma sample generally requires a 2-fold dilution in the 1×Assay buffer. If a sample has a proinsulin level greater than the highest standard, the sample should be diluted with 1×Assay buffer and the assay should be repeated. It is recommended that the users establish their own dilution factors based on the concentration range of their samples.

ASSAY PROCEDURE

It is recommended that all standards and samples be assayed in duplicate.

1. Add 50 µL of 1×Detection antibody solution to each well.
2. Add 50 µL of standard or sample per well. Seal the plate with a plate cover. Incubate at room temperature for 1 hour, shaking the plate at 600 rpm on a horizontal micro-plate shaker.
3. Discard the content and tap the plate on a clean paper towel to remove residual solution in each well. Add 300 µL of 1×Wash buffer to each well and incubate for 1 minute. Discard the 1×Wash buffer and tap the plate on a clean paper towel to remove residual wash buffer. Repeat the wash step for a total 5 washes.
4. Add 100 µL of Substrate solution to each well, incubate at room temperature for 15 minutes. **Protect from light.**
5. Add 100 µL of Stop solution to each well, gently tap the plate frame for a few seconds to ensure thorough mixing.
6. Measure absorbance of each well at 450 nm immediately.

CALCULATION

1. Subtract the absorbance of the blank from that of standards and samples.
2. Generate a standard curve by plotting the absorbance obtained (y-axis) against human proinsulin concentrations (x-axis). The best fit line can be generated with any curve-fitting software by regression analysis. Any curve of 4-parameter or log-log curve fitting can be used for calculation.
3. Determine human proinsulin concentration of samples from standard curve.

Website: www.immunodiagnostics.com.hk

E-mail: info@immunodiagnostics.com.hk(HK) /info@immunodiagnostics.ca(Canada)

Tel: +852 3502 2780 (HK) / +1-437-886-5136 (Canada)


Version 09/25



TYPICAL STANDARD CURVE

The following standard curve is provided for demonstration only. A standard curve should be generated for each set of sample assay.

Human proinsulin (pg/mL)	Absorbance (450 nm)	Blanked Absorbance
0	0.087	0
31.2	0.107	0.020
62.5	0.141	0.054
125	0.223	0.136
250	0.374	0.287
500	0.794	0.707
1000	1.494	1.407
2000	2.599	2.512

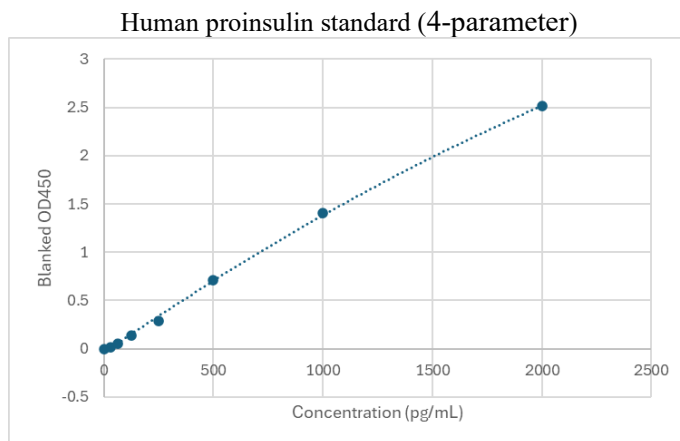

ASSAY CHARACTERISTICS

A. Sensitivity

The lowest level of proinsulin that can be measured by this assay is 31.2 pg/mL.

B. Precision

Intra-assay Precision (Precision within an assay) C.V. <6.5%.

Inter-assay Precision (Precision between assays) C.V. <7.3%.

Website: www.immunodiagnostics.com.hk

E-mail: info@immunodiagnostics.com.hk(HK) /info@immunodiagnostics.ca(Canada)

Tel: +852 3502 2780 (HK) / +1-437-886-5136 (Canada)

Version 09/25



SUMMARY OF ASSAY PROCEDURE

Add 50 μ L of 1 \times Detection antibody solution to each well.



Add 50 μ L of standard or sample to each well.



Incubate at room temperature for 1 hour (600 rpm).



Aspirate and wash each well five times.



Add 100 μ L of Substrate solution to each well.



Incubate at room temperature for 15 minutes.



Add 100 μ L of Stop solution to each well.



Measure absorbance of each well at 450 nm.



Calculation

Website: www.immunodiagnostics.com.hk

E-mail: info@immunodiagnostics.com.hk(HK) / info@immunodiagnostics.ca(Canada)

Tel: +852 3502 2780 (HK) / +1-437-886-5136 (Canada)

Version 09/25